HI96832

Propylene Glycol Refractometer

NSTRUCTION MANUAL





Dear Customer,

Thank you for choosing a Hanna Instruments product.

Please read this instruction manual carefully before using this instrument.

This manual will provide you with the necessary information for correct use of this instrument, as well as a precise idea of its versatility.

If you need additional technical information, do not hesitate to e-mail us at tech@hannainst.com or view our worldwide contact list at www.hannainst.com.

All rights are reserved. Reproduction in whole or in part is prohibited without the written consent of the copyright owner, Hanna Instruments Inc., Woonsocket, Rhode Island, 02895, USA.

Preliminary Examination4
General Description5
Specifications6
Principle of Operation7
Measurement Units8
Functional Description9
Display Elements10
Measurement Guidelines
Calibration Procedure
Measurement Procedure
To Change Measurement Unit14
To Change Temperature Unit14
Making a Standard Solution15
Battery Replacement
Error Messages
Accessories

Remove the instrument from the packing materials and examine carefully to ensure no damage has occurred during shipping. If any damage has occurred, contact your local Hanna Instruments Office.

Each H196832 instrument is supplied with:

- 9 V battery
- Instruction manual
- Plastic pipette
- Quality Certificate

Note: Save all packing material until you are sure that the instrument functions correctly. A defective instrument must be returned in its original packing.

The H196832 Digital Refractometer is a rugged portable, water resistant device that utilizes the measurement of refractive index to determine the % volume and freezing point of propylene glycol based solutions.

Propylene glycol is used in many applications including:

- pharmaceutical solvent
- · humectant food additive
- emulsification agent
- moisturizer in cosmetics, toothpaste, tabacco products, etc.
- hand sanitizer and lotion
- antifreeze

The H196832 device benefits from Hanna Instruments' years of experience as a manufacturer of analytical instruments. The digital refractometer eliminates the uncertainty associated with mechanical refractometers and is easily portable for use in the field.

The HI96832 refractometer is an optical device that is simple and quick to use. Samples are measured after a simple user calibration with distilled or deionized water. Within seconds, the refractive index and temperature are measured and converted into one of two measurement units: % Volume or Freezing Point. The instrument utilizes internationally recognized references for unit conversion and temperature compensation for propylene glycol solutions (e.g. CRC Handbook of Chemistry and Physics, 87th Edition).


The temperature (in ${}^{\circ}\text{C}$ or ${}^{\circ}\text{F}$) is also displayed on the large dual level display along with helpful message codes.

Key features include:

- IP65 waterproof protection
- Automatic Temperature Compensation (ATC)
- Battery operation with Low Power indicator (BEPS)
- Automatically turns off after 3 minutes of non-use

	% Volume: 0.0 to 100.0 %V/V		
Range	Freezing Point: 0.0 to -51.0 °C (32.0 to -59.8 °F)		
	°C (°F): 0.0 to 80.0 °C (32.0 to 176.0 °F)		
	% Volume: 0.1 %		
Resolution	Freezing Point: 0.1 °C (0.1 °F)		
	°C (°F): 0.1 °C (0.1 °F)		
Accuracy	$\%$ Volume: $\pm 0.3~\%$		
	Freezing Point: ± 0.5 °C (± 1.0 °F)		
	°C (°F): ±0.3 °C (±0.5 °F)		
Temperature Compensation	Automatic between 0.0 and 40.0 °C (32.0 - 104.0 °F)		
Measurement Time	Approximately 1.5 seconds		
Minimum Sample Volume	100 μL (cover prism totally)		
Light Source	Yellow LED		
Sample Cell	Stainless Steel ring and flint glass prism		
Case Material	ABS		
Enclosure Rating	IP65		
Battery Type / Life	9V / 5000 readings		
Auto-Off	After 3 minutes of non-use		
Dimensions	192 x 102 x 69 mm (7.6 x 4.1 x 2.7")		
Mass	350 g (12.3 oz.)		


Propylene Glycol determinations are made by measuring the refractive index of solution. Refractive Index is an optical characteristic of a substance and the number of dissolved particles in it. Refractive Index is defined as the ratio of the speed of light in empty space to the speed of light in the substance. A result of this property is that light will "bend", or change direction, when it travels through a substance of different refractive index. This is called refraction. When passing from a material with a higher to lower refractive index, there is a critical angle at which an incoming beam of light can no longer refract, but will instead be reflected off the interface.

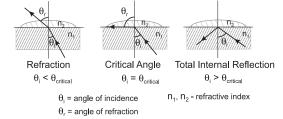

The critical angle can be used to easily calculate the refractive index according to the equation:

$$\sin{(\theta_{\text{critical}})} = n_{2}/n_{1}$$

Where ${\bf n_2}$ is the refractive index of the lower-density medium; ${\bf n_1}$ is the refractive index of the higher-density medium.

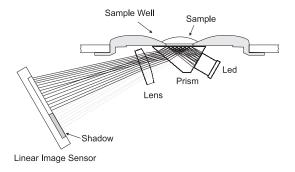


In the H196832 refractometer, light from an LED passes through a prism in contact with the sample. An image sensor determines the critical angle at which the light is no longer refracted through the sample.


Specialized algorithms then apply temperature compensation to the measurement and convert the refractive index to: % Volume or Freezing Point.

Freezing point is displayed as a temperature from 0.0 to -50.0 $^{\circ}\text{C}$ corresponding to 0.0-60.0 % by volume.

When the display shows "LO", the freezing point is at a minimum (below -51 °C).

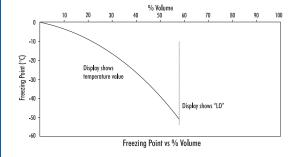


TOP VIEW

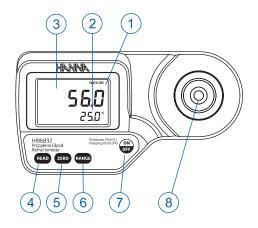


- 1. Liquid Crystal Display (LCD)
- 2. Secondary Display
- 3. Primary Display
- 4. READ Key (User Measurement)
- 5. ZERO Key (User Calibration)
- 6. RANGE Key (User Measurement Unit)
- 7. **ON/OFF**
- 8. Stainless Steel Sample Well and Prism

BOTTOM

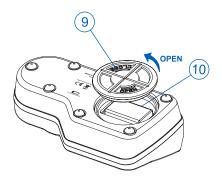

- 9. Battery Cover
- 10. Battery Compartment

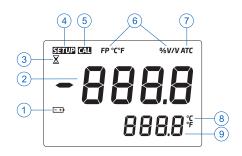

- 1. Battery (blinks when low battery condition detected)
- 2. Primary Display (displays measurement and error messages)
- 3. Measurement in Progress Tag
- 4. SETUP: Factory Calibration Tag
- 5. Calibration Tag
- 6. Measurement Unit
- 7. Automatic Temperature Compensation (blinks when temperature exceeds 0.0-40.0 °C / 32.0-104.0 °F range)
- 8. Temperature Units
- 9. Secondary Display (displays temperature measurements; when blinking, temperature has exceeded operation range: 0.0-80.0 °C/32.0-176.0 °F)

- Handle instrument carefully. Do not drop.
- Do not immerse instrument under water.
- Do not spray water to any part of instrument except the "sample well" located over the prism.
- The instrument is intended to measure propylene glycol solutions. Do not expose instrument or prism to solvents that will damage it. This includes most organic solvents and extremely hot or cold solutions.
- Particulate matter in a sample may scratch the prism. Absorb sample with a soft tissue and rinse sample well with deionized or distilled water between samples.
- Use plastic pipettes to transfer all solutions. Do not use metallic tools such as needles, spoons or tweezers as these will scratch the prism.
- Cover sample well with hand if measuring in direct sun.
- To reduce the effects of evaporation or absorption of water when taking readings over a period of time, the prism and sample well can be covered with plastic wrap.

Calibration should be performed daily, before measurements are made, when the battery has been replaced, between a long series of measurements, or if environmental changes have occurred since the last calibration.

 Press the ON/OFF key, then release. Two instrument test screens will be displayed briefly; all LCD segments followed by the percentage of remaining battery life. The meter will then briefly display an indication of the measurement units set. When LCD displays dashes, the instrument is ready.





2. Using a plastic pipette, fill the sample well with distilled or deionized water. Make sure the prism is completely covered.



Note: If the ZERO sample is subject to intense light such as sunlight or another strong source, cover the sample well with your hand or other shade during the calibration.

3. Press the **ZERO** key. If no error messages appear, your unit is calibrated. (For a description of ERROR MESSAGES see page 17).







Note: The 0.0 screen will remain until a sample is measured or the power is turned off.

 Gently absorb the ZERO water standard with a soft tissue. Use care not to scratch the prism surface. Dry the surface completely. The instrument is ready for sample measurement.



Note: If the instrument is turned off, the calibration will not be lost.

Verify the instrument has been calibrated before taking measurements.

 Wipe off prism surface located at the bottom of the sample well. Make sure the prism and sample well are completely dry.



Using a plastic pipette, drip sample onto the prism surface. Fill the well completely.



Note: If the temperature of the sample differs significantly from the temperature of the instrument, wait approximately 1 minute to allow thermal equilibration.

3. Press the **READ** key. The results are displayed as % Volume or Freezing Point.







Note: The last measurement value will be displayed until the next sample is measured or the instrument is turned off. Temperature will be continuously updated.

The ATC tag blinks and automatic temperature compensation is disabled if the temperature exceeds the 0.0-40.0 $^{\circ}$ C / 32.0-104.0 $^{\circ}$ F range.

- 4. Remove sample from the sample well by absorbing with a soft tissue.
- Using a plastic pipette, rinse prism and sample well with distilled or deionized water. Wipe dry. The instrument is ready for the next sample.



Press the **RANGE** key to select the measurement units. The instrument toggles between the two measurement scales each time the key is pressed and the primary display indicates "**VoL**" % Volume or "FrPt" for Freezing Point. When the instrument displays the screen with 4 dashes, the instrument is ready for measurement.



To change the temperature measurement unit from Celsius to Fahrenheit (or vice versa), follow this procedure.

 Press and hold the ON/OFF key continuously for approximately 8 seconds. The LCD will display the "all segment" screen followed by a screen with the model number on the primary display and the version number on the secondary display. Continue pressing the ON/OFF key.



2. While continuing to hold the **ON/OFF** key, press the **ZERO** key. The temperature unit will change from °C to °F or vice versa.



Note: The selected unit will be used in Freezing Point mode.


To make a Propylene Glycol Solution, follow the procedure below:

- To make an X % propylene glycol solution volumetrically add X mL of high purity propylene glycol (CAS #: 57-55-6; MW 76.09) to a 100 mL Grade A volumetric flask.
- Use distilled or deionized water to bring the flask close to the total volume, mix, allow the solution to return to room temperature.
- Once the solution has returned to room temperature use distilled or deionized water to bring the total volume to 100 mL. Mix solution well before use.

	Propylene Glycol	Total Volume	Expected Freezing Point Value
10 % V	10.00 mL	100.00 mL	-3.4 °C (25.9 °F)
40 % V	40.00 mL	100.00 mL	-21.3 °C (-6.3 °F)

To replace the instrument's battery, follow these steps:

- Make sure the instrument is off.
- Turn instrument upside down and remove the battery cover by turning it counterclockwise.

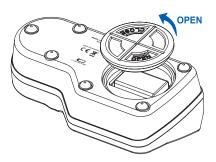

- Extract the battery from its location.
- Replace with a new 9V battery making sure to observe polarity.
- Insert the back battery cover and fasten it by turning clockwise to engage.

Error Code		Description
"Err"	Err 25.0°	General failure. Cycle power to instrument. If error persists, contact your local Hanna Instruments Office.
"LO" primary display	L 0 25.0°	Sample exceeds minimum measurement range.
"HI" primary display	HI 25.0°	Sample exceeds maximum measurement range.
"LO" primary display Cal segment ON	L 0 25.0*	Wrong solution used to zero instrument. Use deionized or distilled water. Press ZERO.
"HI" primary display Cal segment ON	H ;	Wrong solution used to zero instrument. Use deionized or distilled water. Press ZERO.
"t LO" primary display Cal segment ON	E LO 0.0°	Temperature exceeds ATC low limit (10.0 °C) during calibration.
"t HI" primary display Cal segment ON	E HI 40.3*	Temperature exceeds ATC high limit (40.0 °C) during calibration.
"Air"	A, r	Prism surface insufficiently covered.
"ELt"	ELE 25.0°	Too much external light for measurement. Cover sample well with hand.
"nLt"	nLE 25.0°	LED light is not detected. Contact your local Hanna Instruments Office.
Battery segment blinking	56.0	<5% of battery life is remaining.
Temperature values are blinking "0.0°C" or "80.0°C"	5 6.0	Temperature measurement out of range (0.0 to 80.0 °C).
ATC segment blinking	56.0	Outside temperature compensation range (10.0 to 40.0 °C).
SETUP segment blinking	 25.0°	Factory calibration lost. Contact your local Hanna Instruments Office.

Code	Description
HI740157P	Plastic refilling pipette (20 pcs.)
HI740029P	9V battery (10 pcs.)

Warranty

H196832 is warranted for two years against defects in workmanship and materials when used for their intended purpose and maintained according to instructions.

This warranty is limited to repair or replacement free of charge.

Damage due to accidents, misuse, tampering or lack of prescribed maintenance is not covered.

If service is required, contact your local Hanna Instruments Office. If under warranty, report the model number, date of purchase, serial number and the nature of the problem. If the repair is not covered by the warranty, you will be notified of the charges incurred. If the instrument is to be returned to Hanna Instruments, first obtain a Returned Goods Authorization number from the Technical Service department and then send it with shipping costs prepaid. When shipping any instrument, make sure it is properly packed for complete protection.

To validate your warranty, fill out and return the enclosed warranty card within 14 days from the date of purchase.

Hanna Instruments reserves the right to modify the design, construction or appearance of its products without advance notice.

World Headquarters

Hanna Instruments Inc. Highland Industrial Park 584 Park East Drive Woonsocket, RI 02895 USA www.hannainst.com

Local Office

Hanna Instruments USA 270 George Washington Highway Smithfield, RI 02917

Phone: 800.426.6287 Fax: 401.765.7575

e-mail: tech@hannainst.com

